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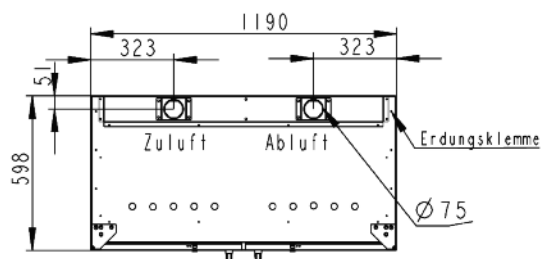
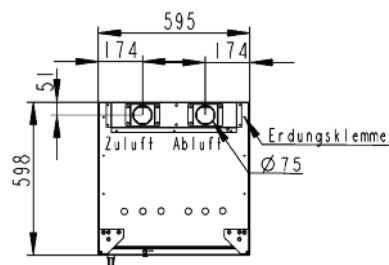
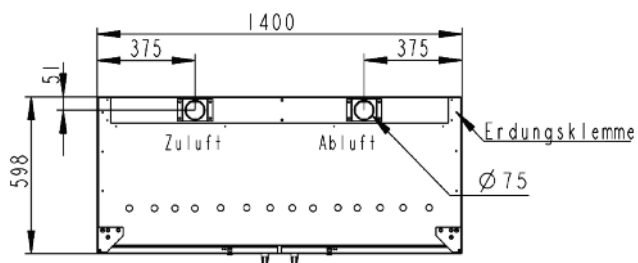
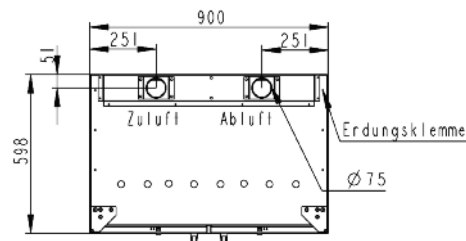
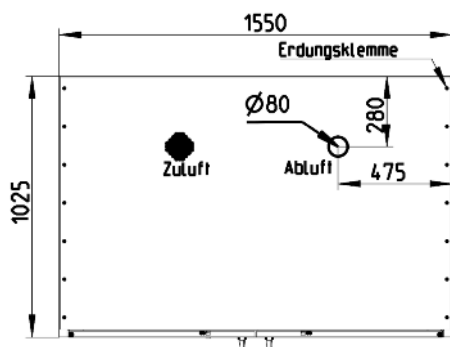
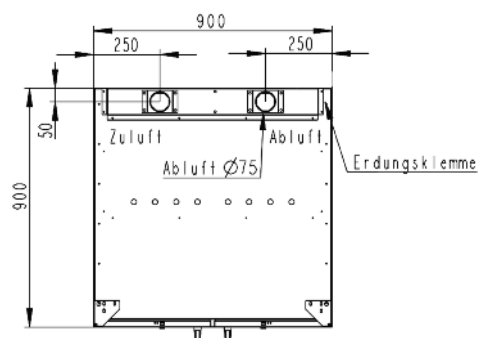
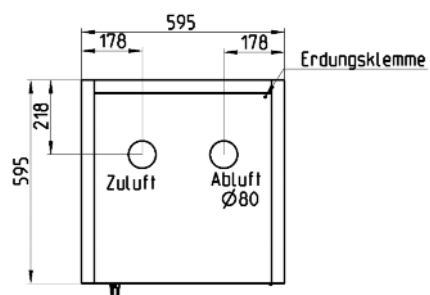
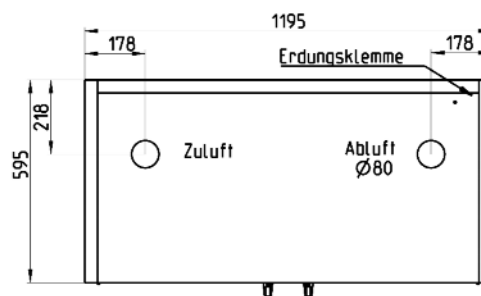
SiZ G30

[illegible]

Technical drawing of a rectangular plate with the following dimensions and features:

- Overall width: 595
- Overall height: 506
- Inner width: 473
- Inner height: 428
- Top-left corner: Zulauf (Inlet)
- Top-right corner: Ablauf (Outlet)
- Top-right corner detail: Entlastungskehle (Relief notch)
- Bottom-left corner: Ø 25 (Hole)
- Bottom edge: 5x Ø 25 (5 holes)
- Center hole: Ø 100

Technical drawing of a rectangular ventilation unit. The overall dimensions are 1400 mm in width and 598 mm in height. The unit features a 'Zuluft' (supply air) inlet on the left and an 'Abluft' (exhaust air) outlet on the right. The distance between the centerlines of these openings is 293 mm. The unit is equipped with 16 circular fans, each with a diameter of $\varnothing 25$ mm, arranged in a single row along the bottom. The distance between the centerlines of these fans is 80 mm. The unit is connected to a power supply via an 'Erddungsklemme' (grounding terminal) on the right side. The distance from the top edge to the centerline of the exhaust air outlet is 156 mm, and the distance from the bottom edge to the centerline of the exhaust air outlet is 428 mm. The unit is shown in a perspective view with a top-down and side view.

Technical drawing of a square ventilation grille. The overall dimensions are 570 mm by 570 mm. The central circular opening has a diameter of $\varnothing 80$ mm. The distance from the center of the opening to the top edge is 196 mm. Labels indicate 'Zuluft' (supply air) on the left, 'Abluft' (exhaust air) on the top, and 'Erdungsklemme' (grounding terminal) on the right. The drawing shows a square frame with mounting points at the corners and midpoints of the sides.

Technical drawing of the front view of a rectangular enclosure. The drawing shows a rectangle with a width of 1100 and a height of 570. On the top edge, there is a circular feature with a diameter of $\varnothing 80$. Above this feature, there is a dimension of 196. Labels with arrows point to the top edge (Zuluft), the circular feature (Abluft), and a terminal (Erduungsklemme).

Technical drawing of a rectangular box. The overall width is 877 and the overall height is 570. A top flange has a width of 195. A circular feature on the top flange has a diameter of $\varnothing 80$. Labels with arrows point to 'Zuluft' (top left), 'Abluft' (top center), and 'Erddungsklemme' (top right).

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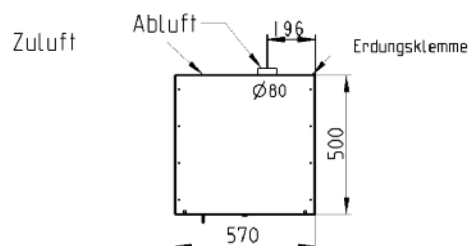
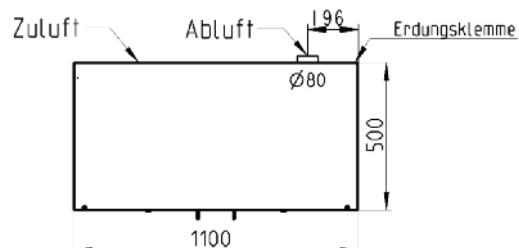
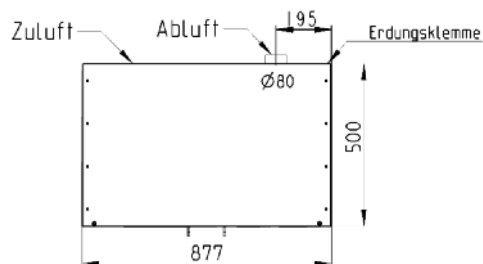
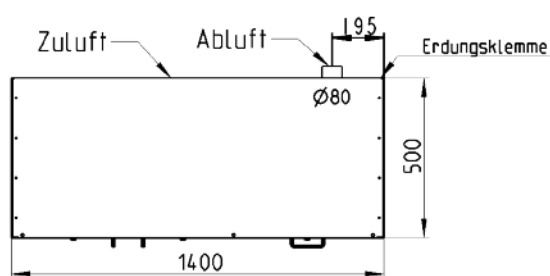
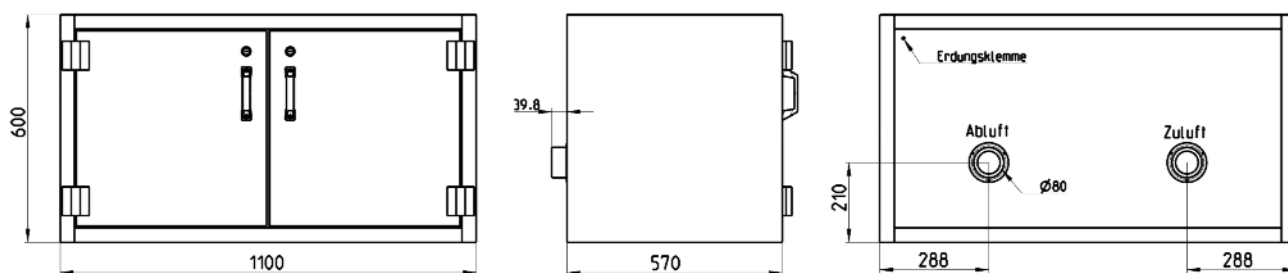
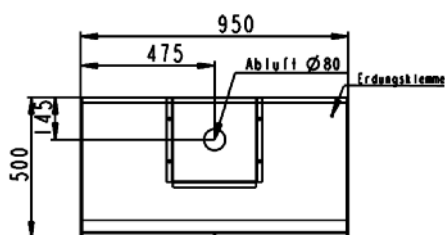
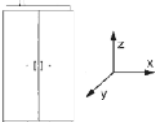
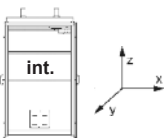
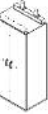
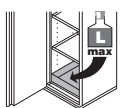
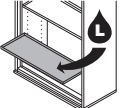
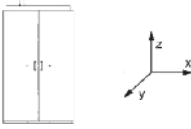
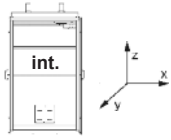
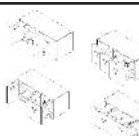
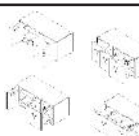
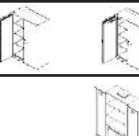
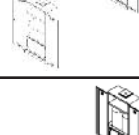
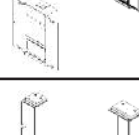
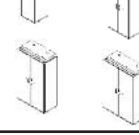
**storeLAB® Safety Storage-Cabinet model 1400 / model 1200 / model 900 /
model 600 / model SIS-FAS 1550 / model SIS-FAS 900
Underbench cabinets model 1400 FT / model 1400 SF / model 1400 SFSF /
model 1400 FTSF /
model 1100 FT / model 1100 SF / model 1100 SFSF / model 1100 FTSF /
model 880 FT / model 880 SF / model 880 FT SF / model 880 SF SF /
model 600 FT / model 600 SF
model 1400-500FT / model 1400-500 SF / model 1400-500 FT SF / model 1400-500SFSF
model 1100-500 FT / model 1100-500 SF / model 1100-500 FT SF /
model 1100-500 SF SF /
model 880-500 FT / model 880-500 SF / model 880-500 FT SF / model 880-500 SF SF /
model 600-500 FT / model 600-500 SF /
Safety cabinets for documents (DKS) model 1200 / model 1400
Chemical cabinets with safety box model 950-SiB 30 / model 950 GL-SiB 30 /
model 950-SiB 60 / model 950 GL-SiB 60 / model 950-2FAS-SiB 30 / model 950-2FAS
GL-SiB 30 / model 950-2FAS-SiB 60 / model 950-2FAS GL-SiB 60**

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to which this declaration relates is in conformity with the relevant provisions of the following stand-
ards or other normative documents.

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DIN EN 14470-1 (2004-07)

Fire safety storage cabinets – Part 1: Safety storage cabinets for
Flammable liquids

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DIN/ISO 3864 Safety colours and safety signs

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Applied nation-
al standards
(germany)

DIN 4844-1 Safety colours and safety signs

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Following the provisions of Directives (Germany)

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and the regulations of the Statutory Accident Insurance Institution

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The operation company is responsible for installing the products in accordance to national regulations
which may apply in respect to the installation, use and also for regular inspections.

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to which this declaration relates is in conformity with the relevant provisions of the following standards or other normative documents.

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EG-Maschinenrichtlinie 2006/42/EG, L'157/65 vom 9.6.2006
EEC-Machinery Directive 2006/42/EEC
EN 292 Sicherheit von Maschinen; Grundbegriffe, allgemeine
Gestaltungsgrundsätze
EN 292 Safety of machinery – Basic concepts, general principles
For design

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